INGLOBAMENTO / MOUNTING

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Generally, the englobing of the sample is necessary for these reasons:

The sample is small or it has an irregular shape therefore the handling is difficult.

It is necessary to examine the edge of the sample for the analysis of the thickness of electrochemical coatings, decarburization, heat treated surface, performing microhardness, etc. ...

Facilitate, under certain conditions, the identification of the sample and archiving.

For englobing the sample, depending on the requirements, it is possible to use the cold resin or the hot resin; for the latter, you have to use a mounting press, which, reaching high temperature, and guaranteeing the necessary pressure, hardens the resin making the metallographic sample.

Hot mounting allows to obtain a planar parallel sample. Hot resins are suitable for samples capable of withstanding high pressure and temperatures ($\sim 175^\circ$)

The cold resin consists in two or more components mixing together (resin + liquid) which are poured into a mold of appropriate size in which the sample is placed. Hardening takes place by chemical reaction with evolution of heat.

Cold resins are suitable for fragile and sensitive samples, which cannot be subjected to high pressure and temperatures.

On the other hand, vacuum cold mounting eliminates any possible air bubbles.

In porous materials, vacuum allows the resin to penetrate into the micro-cavities, compacting the specimen and improving it during the polishing phase.

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You can minimize shrinkage and improve edge retention by cooling the mount to room temperature under pressure before removing it from the mounting press

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Uncured mounts can be caused by too much moisture in the mounting compound.
Make sure to properly close the container between uses.

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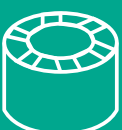
Radial splitting of mounts is often caused by sharp edges on the samples or by samples that are too large for the mold.
Round off sharp corners and move the specimen farther from the edge of the mount.

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Bulging or soft mounts are caused by insufficient cure times. Increase the cure time.

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Unfused or frosted mounting compound is often a sign of insufficient molding temperature or pressure.
Ensure the temperature and pressure settings on the mounting press match the recommendations for the compression compound being using.

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The presence of cracks on the surface of the sample is due to excessive cooling rate.

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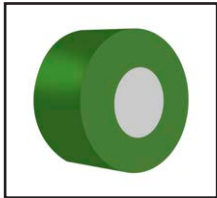
The presence of 'flakes' inside the specimen is due to too low temperatures or the too short duration of the cycle.

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
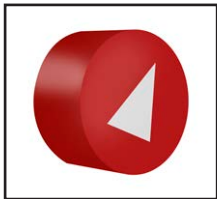
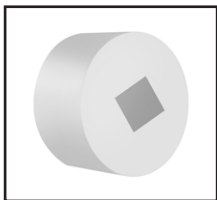
EPOXY RESIN glass fiber filled - EPOGLASS 3

KG	CODE
2	TEC-011-E

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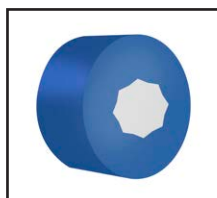DIALLYLPHTHALATE glass fiber filled - GLASS BLU

KG	CODE
1	GG06DAP1000

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Electroconductive thermoplastic resin for electrolytic cleaning and SEM - CONDUCT MOUNT 2


KG	CODE
1	=119906

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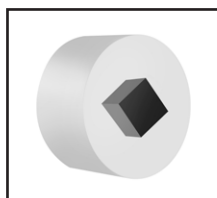
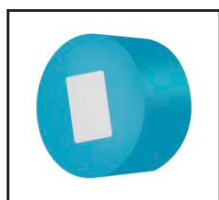
Metacrilato resina TEC-WHITE / TEC-WHITE Methacrilate resin

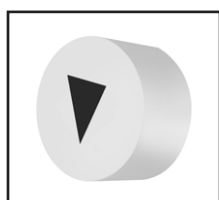
Bianca opaco tricomponente / White opaque three components (1,5Kg powder/1Lt liquid 2/500ml liquid 1)

Pack	Resina / Resin
1,5 kg	=111041

Metacrilato resina TEC-GREEN / TEC-GREEN Methacrilate resin

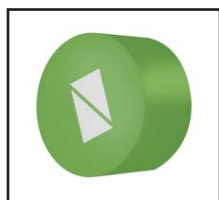
Verde opaco bicomponente (indurimento 5 min.) / Green opaque two components (curing time approx 5 min)

Pack	Resina / Resin
1 Kg	TD-G01P
10 kg	TD-G10P
Pack	Liquido indurente / Hardening liquid
500 ml	TD-G005L
5 lt	TD-G5L

Metacrilato resina TEC- TRASP traslucida / TEC-TRASP Methacrilate resin

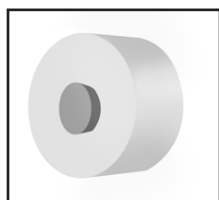Incolore/trasparente bicomponente (indurimento 10 min.) / Colourless/transparent two components (curing time approx 10 min)


Pack	Resina / Resin
1 Kg	TD-T01P
10 Kg	TD-T10P
Pack	Liquido indurente / Hardening liquid
500 ml	TD-T005L
5 lt	TD-T5L

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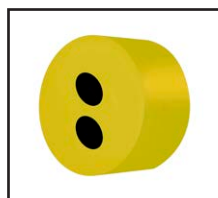Nera opaca bicomponente (indurimento 18 min.) / Black opaque two components (curing time approx 18 min)

Pack	Resina / Resin
1 Kg	TD-K01P
500 ml	TD-K005L

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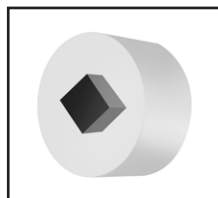
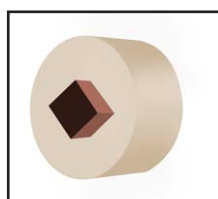
Epoxy Resin low shrinking (curing time approx 15 min) **EPO-FLO**

Pack	Resina / Resin
1 Kg	=111065
250 gr	=111066
1 Kg	=111061
250 gr	=111062

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Multi-pour Vacuum Infiltration Kit

- Struttura a campana
- Rotating plate
- Pompa per il vuoto
- Max capacity
 - 8 silicone mounting cups each Ø 50mm
 - or 8 teflon mounting cups each Ø 40mm.
- incl. 10 mixing cups
- Vacuum pump 230V/50Hz/60W

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REUSABLE MOUNTING MOULDS

	CODICE / CODE	Descrizione / Description
	=113017	Stampini in polypropylene / Polypropylene mould - Ø30 mm, 10 pcs
	=113020	Stampini in polypropylene / Polypropylene mould - Ø40 mm, 10 pcs
	TD-ST25	Stampino in Teflon / Teflon mould Ø 25 mm, 1 pcs
	TD-ST30	Stampino in Teflon / Teflon mould Ø30 mm, 1 pcs
	TD-ST40	Stampino in Teflon / Teflon mould Ø40 mm, 1 pcs
	TD-ST50	Stampino in Teflon / Teflon mould Ø50 mm, 1 pcs
	TD-ST70	Stampino in Teflon / Teflon mould Ø70 mm, 1 pcs
	TD-STR65	Stampino in Teflon / Teflon mould 65x35 mm, 1 pcs
	TD-STR85	Stampino in Teflon / Teflon mould 85x45 mm, 1 pcs
	=113010	Stampo rettangolare in silicone / Rectangular silicon mould 55x30x22 mm, 1 pcs
	=113011	Stampo rettangolare in silicone / Rectangular silicon mould 70x40x22 mm, 1 pcs
	=113012	Stampo rettangolare in silicone / Rectangular silicon mould 100x50x22 mm, 1 pcs
	GG11001	Stampini in silicone / Silicon mould Ø25mm 5 pcs
	GG11002	Stampini in silicone / Silicon mould Ø30mm 5 pcs
	GG11003	Stampini in silicone / Silicon mould Ø1" 1/4 5 pcsL
	GG11004	Stampini in silicone / Silicon mould Ø32mm 5 pcsL
	GG11005	Stampini in silicone / Silicon mould Ø1" 1/2 5 pcs
	GG11006	Stampini in silicone / Silicon mould Ø40mm 5 pcs
	GG11007	Stampini in silicone / Silicon mould Ø50mm 5 pcs

MOLLETTE FERMA CAMPIONI

CLIPS FOR MOUNTING SAMPLES

CODICE / CODE	Descrizione / Description
=113040	Molletta ferma campioni in acciaio inox CLAMP / Specimen support steel clips "CLAMP" for mounting thin samples 100pcs
=113041	Molletta ferma campioni in plastica / Specimen support plastic clips for mounting thin samples 100pcs
=113043	Molletta ferma campioni in plastica nero / Specimen support plastic clips black for mounting thin samples - UNICLIP 100pcs
=113044	Molletta ferma campioni in plastica blu / Specimen support plastic clips blue for mounting thin samples - UNICLIP 100pcs
=113045	Molletta ferma campioni in plastica rosso / Specimen support plastic clips red for mounting thin samples - UNICLIP 100pcs
=113046	Molletta ferma campioni in plastica verde / Specimen support plastic clips green for mounting thin samples - UNICLIP 100pcs
=113047	Molletta ferma campioni in plastica gialla / Specimen support plastic clips yellow for mounting thin samples - UNICLIP 100pcs
=113030	Bicchierini graduati per miscela resine / Graduated mixing cup - 80pcs
=113032	Stick legno per miscela resine / Wooden stirring sticks - 500pcs
FU289200031	Vaschette in polistirene bianco per miscela resine / White polystyrene mixing cups - 90 x 44 x 60 mm 100 pcs
FU289200032	Vaschette in polistirene bianco per miscela resine / White polystyrene mixing cups- 90 x 135 x 25 mm 100 pcs