METALLOGRAPHIC SAMPLE PREPARATION

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The MP 312K Series is designed to offer a highly performing product, and to deliver quality and efficiency at a competitive price.

The innovative design, combined with the base in rigid cast aluminum, and with the cover in shockproof and anti-corrosion material, guarantee its stability while enhancing its functionalities.

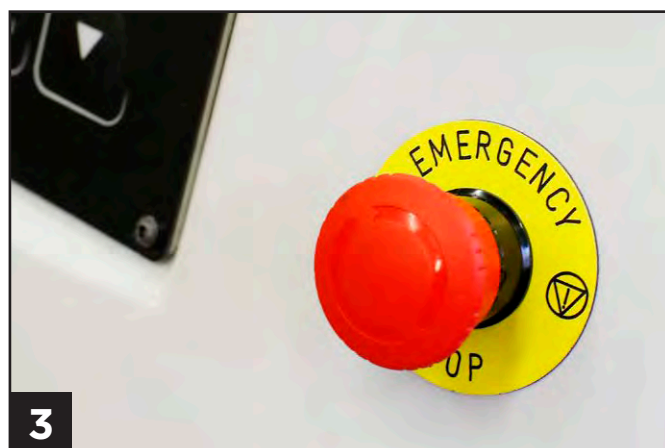
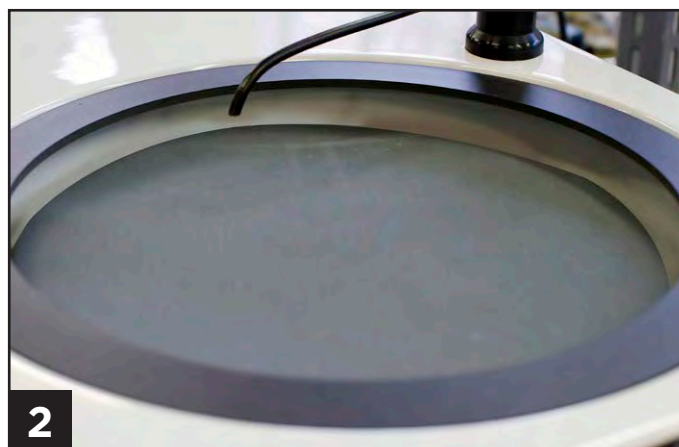
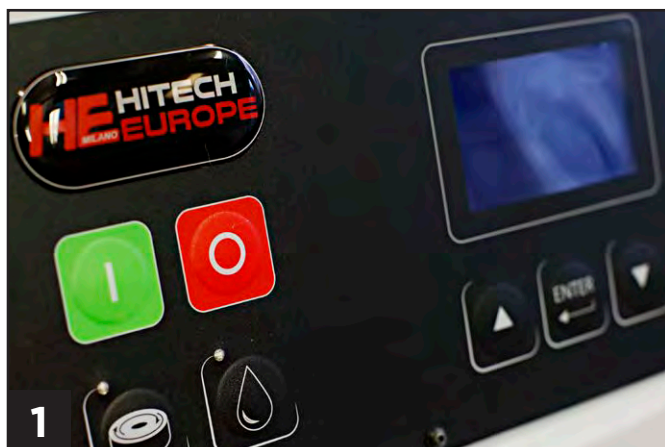
The brushless motor is equipped with an inverter, which optimizes its power, delivering constant torque during the various stages of processing, and maintaining the regularity of rotation.

The electronic board allows to adjust the disc rotational speed (rpm) and direction (clockwise and counterclockwise).

The water flow is managed by a solenoid valve, activated by the electronics that controls its opening and closing, according to the timing set for each grinding and polishing step.

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Display and control panel (1)

Very easy! The front panel consists of a membrane keyboard with the buttons:

- start / stop,
- water tap enabling/disabling
- rotational direction of the working wheel and, of a control panel with digital display for:
- selecting the language
- setting the 'Water' menu
- setting/displaying the working wheel speed rpm and the duration (time) of the cycle
- automatic head settings

Working wheel housing (2)

Suitable for housing working wheels from 200mm up to 300mm diameter, it's been designed with a large capacity, in order to improve and facilitate cleaning operations at the end of the cycle.

Emergency button (3)

Whenever it is necessary to stop the grinding/polishing cycle, by simply pressing the emergency button the working wheel rotation will immediately stop.

Rear panel (4)

The MP 312K Series are preset (in the rear panel) to be connected to peristaltic pump systems for the distribution of fluids, useful for spraying lubricants/diamond suspensions, as well as high density solutions, such as alumina, colloidal silica, etc.

PULITRICI POLISHERS

MP 312KA

- Manual machine
- Single wheel 200/230/250/300 mm
- 0,5Kw Brushless engine with inverter
- Variable speed 0 to 600rpm
- Digital display
- Control panel
- emergency stop button
- Polishing Head and P4K Dispenser preset

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Polishing head and P4K dispenser preset

The new line of 312K series grinder/polishers now also includes the version with the automatic polishing head, which allows you to automatically grind/polish the embedded samples, setting the cycle parameters: air pressure (pneumatic descent of the pistons), duration (timer), water menu. Also have the preset for the digital automatic fluid dispenser P4K.

From the front panel it is possible to set all the parameters to automatically grind/ polish the samples fitted in the sample holder (individual pressure) or in the jig (central pressure):

- number of rpm and direction of rotation of the working wheel and of the head
- Pneumatic pistons descent air pressure
- Cycle duration
- Water ON/OFF

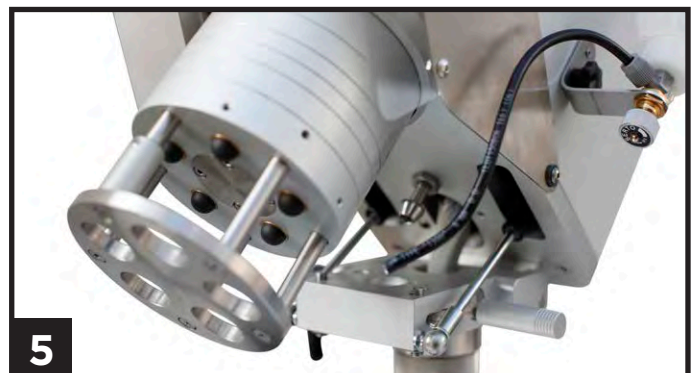
The P4K dispenser, based on peristaltic pumps, is equipped with 4 channels, connected to 4 bottles, allowing you to manage a complete polishing cycle using as needed diamond suspensions, diamond extender (if the suspensions are not diluted) and/or allumina/silica.

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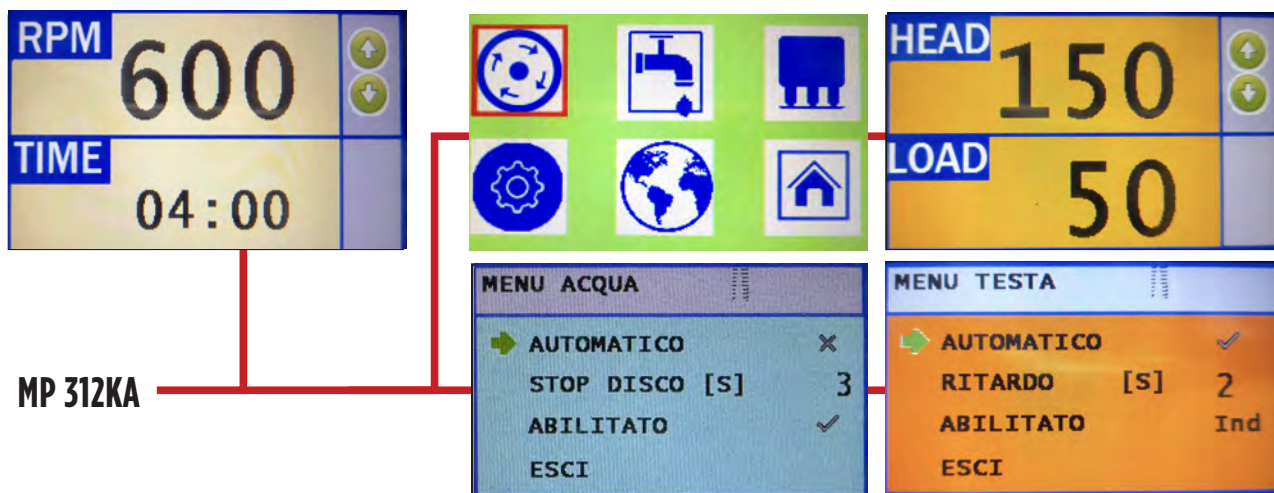
MP 312M - ECONOMIC VERSION

- Manual machine
- Single wheel 200/230/250/300 mm
- 0,3Kw motor
- Variable speed 0 to 600rpm
- front panel:
 - “all-in-one” potentiometer knob (ON/OFF and variable speed setting)
- emergency stop button

ECONOMIC

PULITRICI POLISHERS

MP 312KA

| | |
|--------------------------------------|-------------------------|
| Wheel Ø | From 200 up to 300 mm |
| Brushless Engine | 500W with inverter |
| Engine speed | 0-600 rpm. |
| Head Preset | Yes |
| Dispenser Preset | Yes |
| Rotation | CW - CCW |
| Cycletime | Timed Cycle |
| Power supply | 230V /110V |
| Hydraulic system | Connect to water source |
| Dimensions | 650 x 480 x 240 mm |
| Weight | 38 KG |
| Instruction manual CE certificate | Yes |

POLISHING HEAD

| | |
|-----------|-----------------------|
| IP3-UP | Sample holder 3 holes |
| IP5-UP | Sample holder 5 holes |
| IP5-COMBO | Sample holder 5 holes |

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1 - Paper clamp

2 - Magnetic Base

3 - Working wheel aluminium alloy anticorrosive anodized

4 - Metal Plate

MP 312K SERIES

| MP 312K | | MP 312M | |
|--------------------------------------|-------------------------|--------------------------------------|-------------------------|
| Wheel Ø | From 200 up to 300 mm | Wheel Ø | From 200 up to 300 mm |
| Brushless Engine | 300W with inverter | Engine | 300W with inverter |
| Engine speed | 0-600 rpm. | Engine speed | 0-600 rpm. |
| Rotation | CW - CCW | Rotation | CW - CCW |
| Cycletime | Timed Cycle | Power supply | 230V /110V |
| Power supply | 230V /110V | Hydraulic system | Connect to water source |
| Hydraulic system | Connect to water source | Dimensions | 650 x 480 x 240 mm |
| Dimensions | 650 x 480 x 240 mm | Weight | 38 KG |
| Weight | 38 KG | Instruction manual CE certificate | Yes |
| Instruction manual CE certificate | Yes | | |

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