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The **AP312KA** is a latest generation automatic polishing machine, designed to produce high quality metallographic specimens, at a competitive price.

The innovative design, combined with the base in rigid cast aluminum, and with the cover in shockproof and anti-corrosion material, guarantees its stability while enhancing its functionalities.

The automatic head for specimen housing is extremely versatile and robust and is completed by a safety guard.

The motor installed on this line of polishers is a brushless type, equipped with an inverter and geared motor, to have great torque and constant speed.

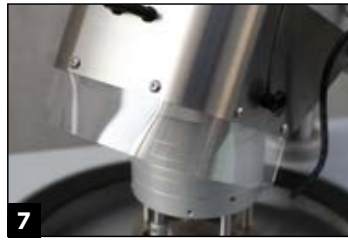
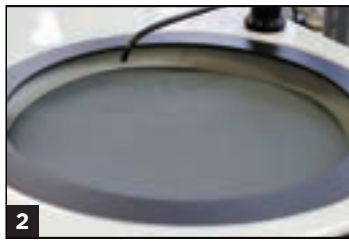
The pneumatic pressure on the specimens is adjusted in a sensitive and precise manner, through a solenoid valve controlled by an electronic board.

The electronic board allows to adjust the disc rotational speed (rpm) and direction (clockwise and counterclockwise).

The water flow is managed by a solenoid valve, activated by the electronics that controls its opening and closing, according to the timing set for each grinding and polishing step.

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Display and control panel (1)

Very easy! The front panel consists of:

- membrane keyboard with the buttons: start / stop, water tap enabling/disabling; rotational direction of the working wheel
- control panel with digital display for: selecting the language; setting the "Water" menu; setting/displaying the working wheel rpm and the duration (time) of the cycle

Working wheel housing (2)

Suitable for housing working wheels from 200mm up to 300mm diameter, it's been designed with a large capacity, in order to improve and facilitate cleaning operations at the end of the cycle.

Emergency button (3)

Whenever it is necessary to stop the grinding/ polishing cycle, by simply pressing the emergency button the working wheel rotation will immediately stop.

Rear panel and preset for Automatic Dispenser (4)

The AP312KA Series are preset (in the rear panel) to be connected to peristaltic pump systems for the distribution of fluids, useful for spraying lubricants/diamond suspensions, as well as high density solutions, such as alumina, colloidal silica, etc.

Head Position Locking Lever (5)

Allows to lock the head in the desired position.

Height adjustment of the head column (6)

The distance between the sample holder and the working wheel can be adjusted by changing the height of the head column using its special flange, depending on the thickness (higher or lower) of the accessories and consumables used during the grinding cycles and polishing.

Safety protection (7)

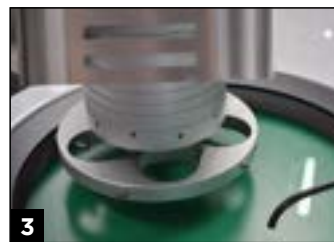
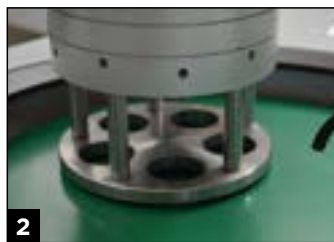
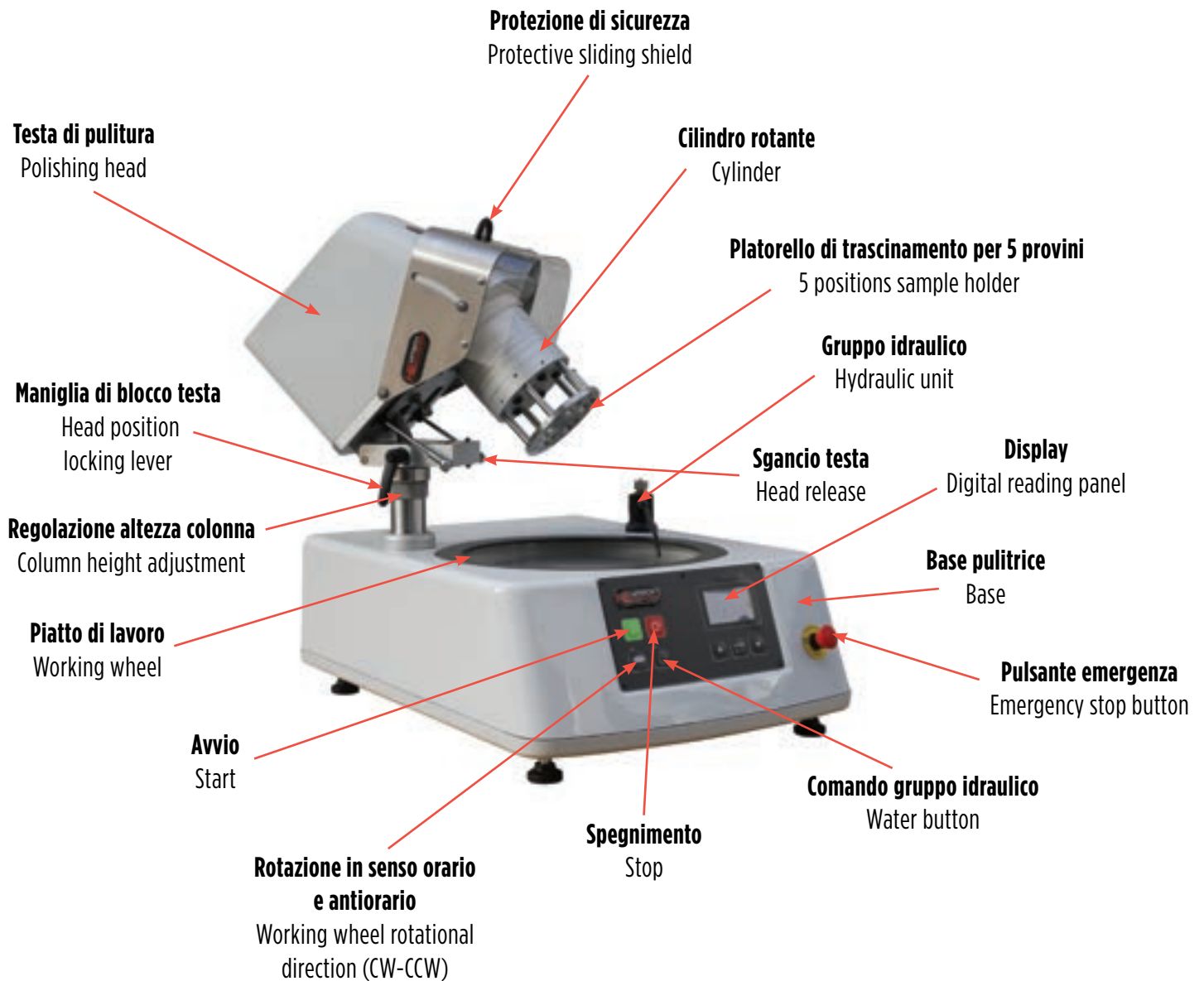
The steel barrier does not allow the operator to bring his hands close to the sample holder while in motion. When the protective shield is shifted down, the safety system gives consent to the rotation of the cylinder or the sample holder, whilst with the shield shifted up the device cannot be used.

Extender manual dispenser (8)

The basic configuration of the machine includes the extender manual dispenser.

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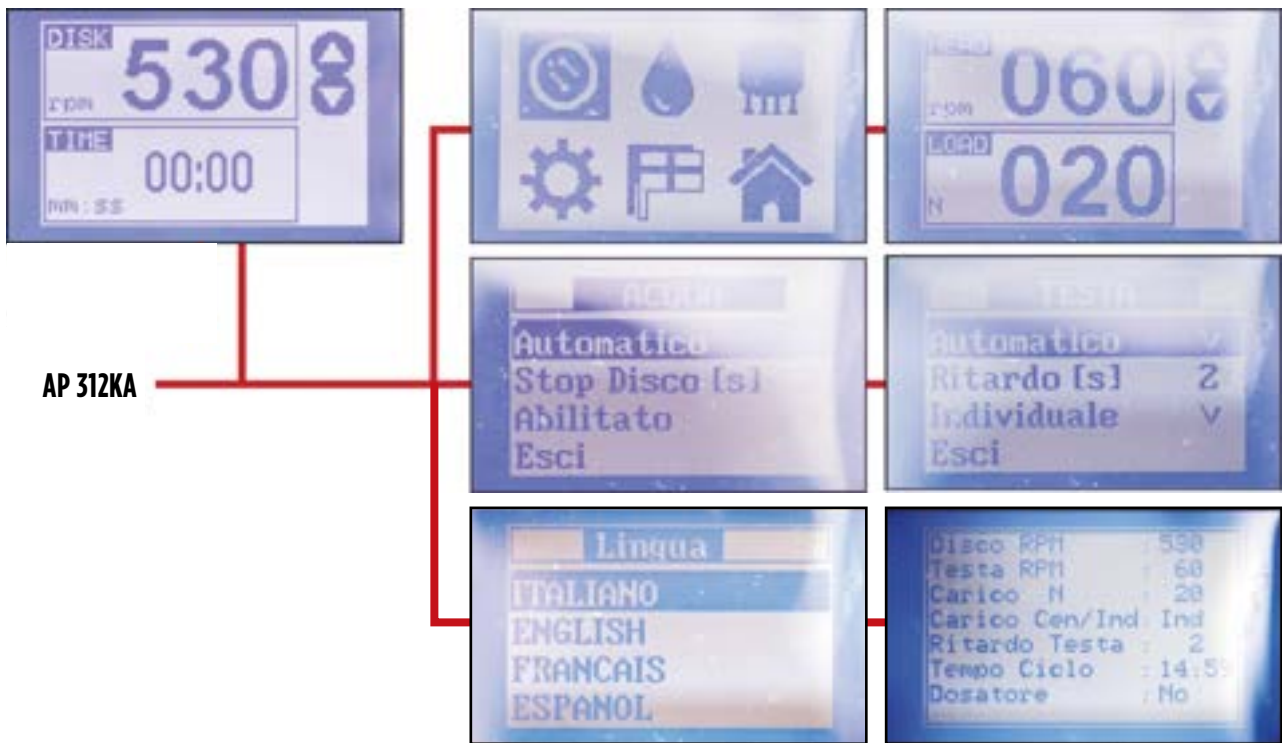
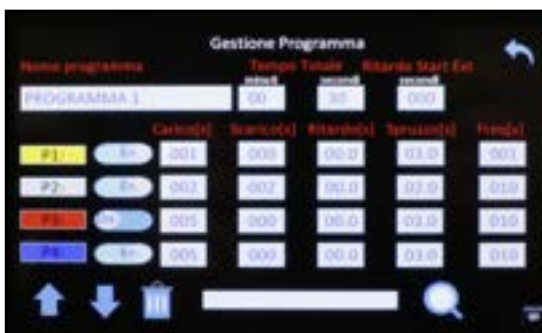
P4K Dispenser, complete with n°4 peristaltic pumps (no. 3 for diluted diamond suspensions and no. 1 also suitable for either colloidal silica or alumina). Touch screen control panel to set cycle parameters and store programs. The P4K can either be connected to the polisher or work separately.

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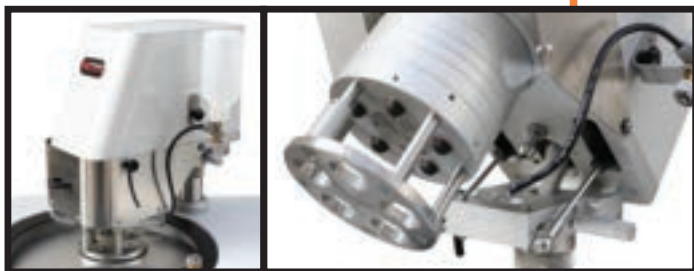
Polishing head and P4 dispenser preset

The new line of 312K series grinder/polishers now also includes the version with automatic polishing head, which allows you to automatically grind/polish the embedded samples, setting the cycle parameters: air pressure (pneumatic descent of the pistons), duration (timer), Water menu. Also have the preset for the digital automatic fluid dispenser P4K.

From the front panel it is possible to set all the cycle parameters to automatically grind / polish the samples fitted in the sample holder (individual pressure) or in the jig (central pressure):

- number of rpm and direction of rotation of the working wheel and of the head
- pneumatic pistons descent air pressure
- cycle duration
- water ON/OFF

The P4K dispenser has 4 peristaltic pumps and 4 channels connected to the 4 bottles, allowing you to manage a complete polishing cycle and use, depending on your needs, diamond suspensions, diamond extender (if the suspensions are not already diluted) and/or alumina/silica.

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AP312KA Polishers – Technical features

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The basic machine does not include working wheel, sample holder, jig/levelling base, sample spacers and paper clamps, to be chosen by user according to the needed sizes.

Sample holders and jigs can be customized.

HE HITECH
MILANO **EUROPE**

METALLOGRAPHIC SAMPLE PREPARATION

tecmet 2000 srl via Monferrato, 5 -20094 Corsico (MI) - Italy
tel+39 02 486 00 385 -tel. +39 02 3663 8452
E mail: tecmet@tecmet2000.it - www.tecmet2000.it