**HE** HITECH  
MILANO **EUROPE**

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The TSP312KA is the latest generation automatic polishing machine, designed to produce high quality metallographic specimens, at a competitive price.

The innovative design, combined with the base in rigid cast aluminum and shockproof, anti-corrosion material, guarantees its stability while enhancing its functionalities.

The automatic head for specimen housing is extremely versatile and equipped with a safety guard.

This line of polishers features a brushless motor, with inverter and geared motor, to have great torque and constant speed.

The pneumatic pressure on the specimens is adjusted in a sensitive and precise manner, through a solenoid valve controlled by an electronic board.

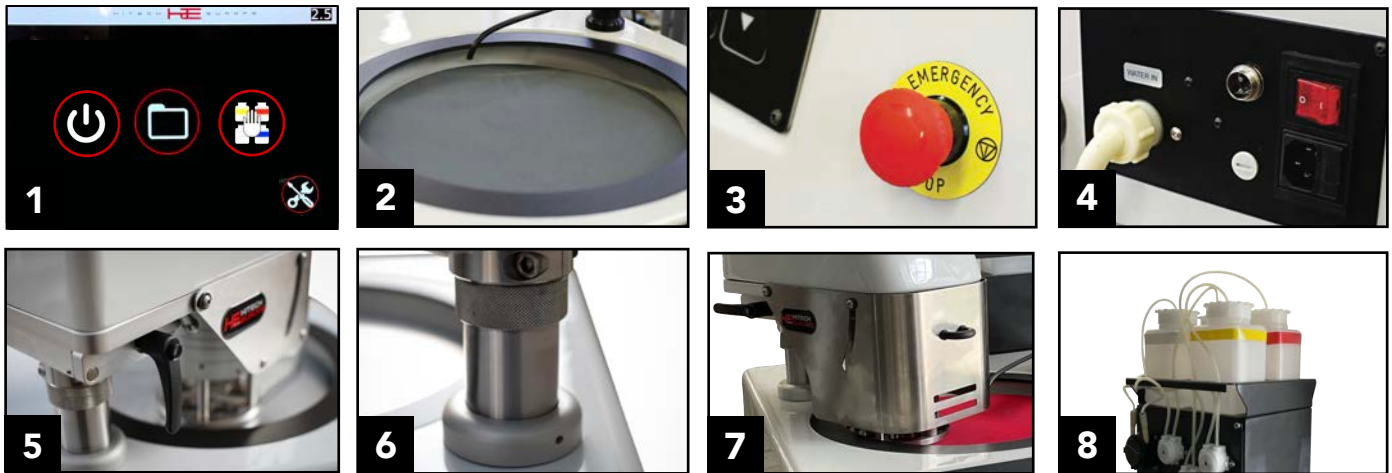
The electronic board allows to adjust the disc rotational speed (rpm) and direction (clockwise and counter-clockwise).

The water flow is managed by a solenoid valve, activated by the electronics that control its opening and closing, according to the timing set for each grinding and polishing step.

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### Touch-screen monitor (1)

Easy-to-use and intuitive touch-screen monitor to adjust all head, plate and pump settings.

### Working wheel housing (2)

Suitable for housing working wheels from 200 mm up to 300 mm diameter, it's been designed with a large capacity, in order to improve and facilitate cleaning operations at the end of the cycle.

### Emergency button (3)

Whenever it is necessary to stop the grinding/polishing cycle, the emergency button will immediately stop the wheel rotation.

### Back control panel (4)

The control panel features: water system connection, air system connection, power supply connection and ON/OFF switch.

### Head position locking lever (5)

Allows to lock the head in the desired position.

### Height adjustment of the head column (6)

The distance between the sample holder and the working wheel can be adjusted by changing the height of the head column using its special flange, depending on the thickness (higher or lower) of the accessories and consumables used during grinding cycles and polishing.

### Safety protection (7)

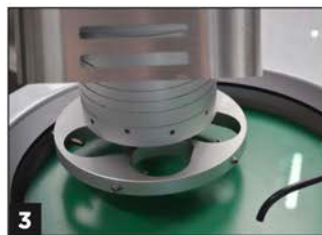
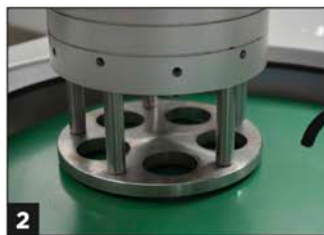
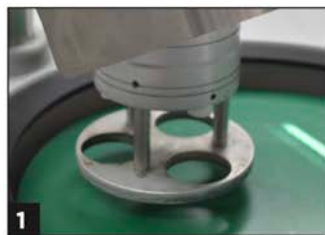
The steel barrier does not allow the operator to bring his hands close to the sample holder while in motion. When the protective shield is shifted down, the safety system gives consent to the rotation of the cylinder or the sample holder, whilst with the shield shifted up the device cannot be used.

### Automatic fluid dispenser (8)

The TSP312KA series features a peristaltic pump system for the distribution of fluids (e.g. lubricants/diamond suspensions) which can be managed through the polisher's touch-screen monitor.

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JIG FOR 4 RECTANGULAR SAMPLES (4)

# TSP 312KA SERIES

## FUNZIONI TOUCH-SCREEN / TOUCH-SCREEN FUNCTIONS

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